

Comprehensive Mycotoxin Analysis: Adding *Alternaria* Toxins and Ergot Alkaloids to Multi-Mycotoxin Analysis in Various Food Commodities

Shun-Hsin Liang*

Restek Corporation, 110 Benner Circle, Bellefonte, PA 16823, USA

ABSTRACT

The ubiquitous contamination of food products with *Alternaria* Toxins (ATs) and Ergot Alkaloids (EAs) often goes unnoticed due to the lack of clear assessments of their human toxicity. To address this issue, a fast and Liquid Chromatography-Mass Spectrometry (LC-MS) friendly method was developed for the accurate and robust analysis of a total of 37 mycotoxins, including ATs and EAs, in various food commodities. This analytical workflow was established through a thorough examination of different types of LC bonding chemistries, sample extraction conditions, and calibration methodologies to ensure accuracy and precision in the measurement.

Keywords: Mycotoxins; *Alternaria* toxins; Ergot alkaloids; *Aspergillus*

DESCRIPTION

Mycotoxins are secondary metabolites produced by several fungal species of the genus *Aspergillus* (aflatoxins and ochratoxins), *Claviceps* (Ergot Alkaloids (EAs)), *Fusarium* (beauvericin, fumonisins, trichothecenes, and zearalenone), *Penicillium* (citrinin, patulin), and *Alternaria* (alternariol, tentoxin, and tenuazonic acid), among others. While ATs and EAs are not currently regulated in the United States, the European Union has established the regulated levels for EAs and is in the process of discussing the maximum permitted levels for ATs [1-3]. EAs contamination is common in cereal crops and wild grasses [4], whereas ATs contamination is prevalent in fruits, vegetables, and cereals [5]. Their co-occurrence in grain-based foods highlights the need for developing an analytical method for the simultaneous analysis of ATs and EAs in such food and feed matrices. However, due to their distinct chemical properties, developing a comprehensive method for simultaneous analysis of multi-class mycotoxins, including emerging ATs and EAs, has been a significant challenge. For liquid chromatography, specific analytical methods using high pH conditions are usually required for the analysis of ATs and EAs to either gain acceptable peak shapes or enable the chromatographic separation of epimers [6-9]. These practices, however, have hindered the simultaneous analysis of these two groups of compounds with other major regulated mycotoxins.

Among more than 70 identified ATs, the five most important ones are altenuene, alternariol, alternariol monomethylether, tentoxin, and tenuazonic acid [10]. Of more than 40 known EAs, the European Union lists six EAs, including ergocornine, ergocristine, ergocryptine, ergometrine, ergosine, and ergotamine, along with their isomeric-inine epimers, as regulated EAs. The guidance is to quantify these 12 EAs separately and sum them up to a total EAs level [11]. Therefore, it is necessary to chromatographically resolve all 12 epimeric EAs for unambiguous and accurate quantification. This study aimed to develop a simplified and unique method for analyzing ATs and EAs together with other major regulated mycotoxins. The method's suitability was evaluated in terms of linearity, specificity, accuracy, precision, and adaptability.

The chromatographic performance was evaluated on Biphenyl, FluoroPhenyl, and C18 stationary phases bonded to superficially porous particles. It was observed that the baseline separation of six EAs and their epimers could only be accomplished on the Biphenyl column under regular acidic conditions.

Similar to other types of multi-compound analysis in food by LC-MS/MS, accurate quantification of multi-mycotoxins in various food commodities requires mitigating the matrix effect. The most direct approach to correct matrix effects is to use stable isotopes as internal standards for quantitative standard calibration. This approach, however, may not be applicable to

Correspondence to: Shun-Hsin Liang, Restek Corporation, 110 Benner Circle, Bellefonte, PA 16823, USA, E-mail: Shun-Hsin.Liang@restek.com

Received: 26-May-2023, Manuscript No. JCT-23-24410; **Editor assigned:** 29-May-2023, PreQC No. JCT-23-24410 (PQ); **Reviewed:** 12-Jun-2023, QC No. JCT-23-24410; **Revised:** 19-Jun-2023, Manuscript No. JCT-23-24410 (R); **Published:** 26-Jun-2023, DOI: 10.35248/2161-0495.23.13.533.

Citation: Liang SH (2023) Comprehensive Mycotoxin Analysis: Adding *Alternaria* Toxins and Ergot Alkaloids to Multi-Mycotoxin Analysis in Various Food Commodities. J Clin Toxicol. 13:533.

Copyright: © 2023 Liang SH. This is an open-access article distributed under the terms of the Creative Commons Attribution License, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited.

most multi-mycotoxin analyses due to the lack of isotopically-labeled compounds for certain analytes. Therefore, the matrix-matched external standard calibration was implemented in this study to equally compensate for the matrix effect of calibration standards and sample solutions for accurate quantification.

Four different food commodities (baby wheat cereal, peanut, tomato puree, and blended flour) were chosen for method validation to demonstrate the applicability of this analytical method across a wide range of food types. Sample extraction was performed using a formic acid-acidified 80:20 Acetonitrile:Water solution followed by extract dry-down and reconstitution in a 50:50 Water:Methanol solution for analysis on a Biphenyl LC column. Chromatographic analysis was performed under LC-MS friendly conditions with a short 11-minute cycle time. The chromatographic and mass spectrometry (Waters Xevo TQ-S) conditions are provided in Table 1, and the resulting chromatogram is shown in Figure 1. Method accuracy and precision were evaluated with samples fortified at 5, 50, and 200 µg/kg. Accurate quantification was achieved using matrix-matched calibration standards at the range of 0.4 to 400 µg/kg. The recoveries of all mycotoxins (except citrinin) in fortified samples were from 70% to 120%, and the relative standard deviation was less than 20% (Table 2). More detailed sample preparation procedures and results can be found in the study of Liang, et al. [12].

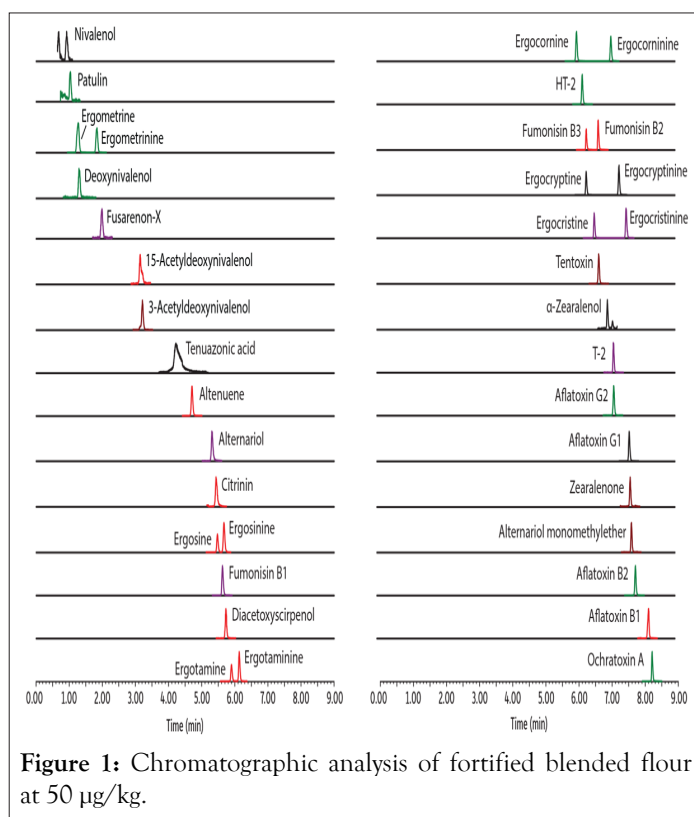


Figure 1: Chromatographic analysis of fortified blended flour at 50 µg/kg.

Chromatographic and mass spectrometry conditions

Analytical column	Raptor Biphenyl 2.7 µm, 100 mm x 2.1 mm	
Guard column	Raptor Biphenyl EXP guard column cartridge 2.7 µm, 5 mm x 2.1 mm	
Mobile phase A	0.05% formic acid in water	
Mobile phase B	0.05% formic acid in methanol	
Gradient	Time (min)	%B
	0	25
	5	50
	9	100
	9.01	25
	11	25
Flow rate	0.4 mL/min	
Injection volume	5 µL	
Column temperature	60°C	
Capillary voltage	0.7 kV	
Gas flow	1000 (L/Hr) Desolvation; 150 (L/Hr) Cone; 7.0 (bar) Nebuliser	
Desolvation temperature	500°C	

Table 1: The chromatographic and mass spectrometry conditions.

Average Recovery (RSD, %)

	Baby wheat cereal			Peanut			Tomato puree			Blended flour		
	Concentration, 5 µg/kg	50	200	5	50	200	5	50	200	5	50	200
Aflatoxin B1	105 (4.8)	100 (3.0)	79.8 (2.6)	98.2 (6.4)	97.0 (5.2)	89.0 (5.7)	92.7 (3.8)	97.6 (5.2)	103 (3.0)	101 (2.8)	95.5 (1.3)	89.0 (1.5)
Aflatoxin B2	110 (1.4)	109 (2.8)	106 (2.3)	102 (5.8)	99.3 (4.7)	91.3 (2.9)	91.7 (4.2)	93.3 (1.9)	94.7 (1.4)	100 (2.3)	101 (1.9)	88.7 (1.3)
Aflatoxin G1	105 (6.1)	107 (1.7)	102 (2.1)	98.2 (4.2)	97.3 (3.2)	91.2 (4.1)	91.3 (1.9)	92.2 (3.6)	93.3 (2.5)	99.3 (1.7)	100 (1.6)	93.6 (2.2)
Aflatoxin G2	108 (3.0)	109 (1.3)	104 (2.2)	104 (5.3)	102 (3.8)	93.5 (1.9)	86.8 (8.3)	96.4 (2.5)	98.5 (2.5)	98.7 (3.1)	102 (2.6)	94.5 (2.0)
Ochratoxin A	109 (1.8)	108 (2.1)	94.5 (1.5)	102 (1.9)	101 (1.1)	97.7 (0.9)	90.9 (3.5)	93.8 (3.3)	101 (5.9)	98.1 (1.6)	98.2 (1.3)	82.8 (1.7)
3- + 15-Acetyldeoxy-nivalenol	104 (6.3)	108 (1.8)	104 (3.3)	101 (6.5)	95.9 (5.8)	91.0 (4.4)	91.9 (4.3)	98.1 (2.7)	95.0 (1.8)	98.4 (5.2)	101 (2.9)	100 (0.9)
Deoxynivalenol	112 (4.0)	102 (2.6)	95.7 (1.3)	98.1 (3.5)	93.7 (4.8)	88.2 (3.4)	-	90.3 (6.4)	94.5 (2.6)	102 (3.5)	97.5 (2.6)	96.9 (1.8)
Diacetoxy-scirpenol	105 (4.0)	107 (1.5)	103 (1.2)	93.2 (4.3)	95.4 (3.9)	93.8 (5.0)	90.9 (3.8)	94.5 (4.7)	94.0 (1.9)	98.1 (6.3)	101 (3.1)	98.7 (1.8)
Fumonisin B1	94.3 (4.6)	94.0 (2.8)	92.3 (2.6)	67.2 (3.1)	88.2 (4.5)	87.8 (6.6)	91.8 (3.6)	91.5 (1.9)	91.9 (0.7)	100 (3.2)	99.6 (1.7)	96.1 (1.2)
Fumonisin B2	93.3 (4.1)	95.1 (4.8)	90.3 (2.9)	95.4 (4.7)	92.5 (2.3)	88.8 (3.9)	89.9 (4.1)	92.9 (2.3)	92.4 (0.8)	104 (2.7)	99.6 (1.4)	94.4 (1.6)
Fumonisin B3	91.8 (4.9)	94.6 (4.9)	91.6 (3.1)	90.6 (2.7)	90.1 (3.8)	87.7 (4.7)	91.1 (3.6)	93.1 (1.8)	91.9 (0.9)	104 (2.2)	99.9 (1.4)	95.9 (1.2)
Fusarenon-X	99.0 (3.9)	100 (2.9)	103 (2.8)	86.9 (7.0)	90.3 (11.0)	88.3 (10.1)	-	92.0 (6.8)	94.3 (1.9)	101 (3.8)	100 (3.7)	98.3 (1.6)
HT-2	110 (2.4)	111 (1.4)	108 (1.1)	100 (2.7)	100 (2.0)	94.3 (3.0)	96.8 (3.1)	96.1 (2.1)	99.0 (1.4)	101 (1.6)	103 (2.2)	98.3 (1.3)
Nivalenol	-	-	-	-	98.3 (6.2)	89.0 (3.6)	-	92.5 (4.5)	93.7 (5.0)	-	95.5 (4.7)	92.9 (2.3)
T-2	111 (2.1)	110 (1.8)	108 (2.8)	99.1 (2.7)	101 (1.7)	95.9 (2.1)	92.0 (6.3)	94.7 (1.3)	98.6 (1.5)	102 (1.3)	103 (1.3)	96.9 (1.3)
α-Zearalenol	100 (4.9)	102 (5.2)	90.1 (5.8)	89.2 (8.1)	93.6 (5.5)	94.7 (3.4)	97.7 (3.2)	88.9 (4.2)	90.0 (3.4)	96.9 (3.7)	99.0 (3.6)	95.0 (3.3)

Zearalenone	110 (6.7)	110 (3.0)	105 (3.7)	98.3 (7.3)	97.4 (2.8)	91.3 (1.5)	95.0 (4.5)	93.6 (2.2)	95.7 (2.0)	101 (3.8)	102 (2.1)	92.3 (1.4)
Citrinin	26.1 (9.2)	26.6 (3.1)	30.1 (3.8)	24.1 (8.7)	25.1 (1.9)	25.8 (3.5)	71.9 (4.7)	76.4 (1.6)	77.1 (1.7)	32.3 (3.5)	32.2 (6.3)	35.8 (4.5)
Patulin	106 (4.6)	95.6 (5.6)	89.2 (5.1)	88.8 (12.0)	83.6 (9.0)	86.0 (7.2)	-	98.9 (3.6)	103 (4.5)	93.6 (4.4)	86.1 (3.1)	92.2 (2.9)
Alternariol	108 (4.1)	108 (1.6)	104 (1.0)	94.2 (3.4)	95.4 (2.4)	96.2 (2.7)	89.3 (4.6)	91.8 (2.5)	91.4 (1.3)	98.4 (2.3)	101 (2.5)	96.3 (3.2)
Alternariol monomethyl ether	108 (4.1)	109 (2.2)	99.3 (2.7)	93.5 (3.3)	93.5 (3.7)	89.8 (2.4)	91.3 (6.6)	88.7 (5.1)	93.9 (3.9)	104 (2.9)	101 (1.7)	93.7 (1.9)
Altenuene	110 (2.1)	109 (2.1)	105 (2.1)	99.6 (2.0)	99.5 (1.2)	95.4 (1.2)	98.4 (3.4)	92.4 (2.1)	92.8 (1.8)	101 (2.9)	101 (3.1)	98.2 (0.5)
Tentoxin	111 (3.6)	109 (2.5)	103 (1.4)	104 (2.9)	101 (1.1)	95.3 (1.4)	92.5 (6.2)	94.2 (2.2)	95.8 (1.4)	104 (4.2)	105 (2.1)	98.2 (1.9)
Tenuazonic acid	-	85.8 (1.7)	87.4 (6.3)	92.5 (4.7)	91.0 (2.1)	88.5 (2.4)	-	89.3 (4.1)	88.5 (2.0)	-	92.5 (8.8)	90.0 (9.5)
Ergocornine	109 (1.5)	109 (1.4)	102 (1.3)	93.8 (3.5)	93.2 (4.4)	91.2 (3.3)	91.5 (3.0)	93.1 (1.9)	92.9 (0.6)	102 (2.5)	101 (1.9)	97.6 (1.7)
Ergocorninine	109 (3.0)	109 (2.0)	101 (1.9)	105 (3.0)	104 (2.4)	99.5 (3.1)	89.9 (3.8)	92.3 (2.2)	92.5 (3.1)	101 (2.5)	102 (2.6)	95.7 (2.4)
Ergocristine	108 (3.1)	108 (2.9)	101 (4.4)	92.1 (3.8)	91.7 (5.1)	92.0 (2.2)	91.3 (2.9)	94.2 (2.0)	94.3 (0.8)	101 (1.7)	99.8 (2.0)	96.7 (1.8)
Ergocristinine	106 (3.5)	105 (1.4)	101 (0.8)	102 (4.8)	104 (4.3)	102 (4.6)	91.6 (5.9)	94.4 (1.8)	95.6 (2.7)	102 (2.9)	102 (3.0)	99.3 (4.5)
Ergocryptine	107 (2.0)	109 (1.9)	104 (3.4)	95.0 (3.0)	94.7 (4.1)	92.1 (1.7)	90.1 (3.0)	93.5 (2.2)	93.2 (0.7)	99.5 (2.7)	99.9 (1.2)	97.4 (1.4)
Ergocryptinine	106 (1.7)	108 (2.0)	101 (1.1)	103 (5.3)	105 (4.0)	101 (4.2)	91.1 (4.3)	95.1 (1.5)	98.1 (1.6)	101 (2.0)	101 (1.8)	95.4 (1.9)
Ergometrine	92.8 (7.3)	90.0 (4.2)	88.3 (3.6)	101 (2.3)	96.2 (2.6)	86.7 (1.9)	90.7 (3.6)	88.9 (6.1)	87.6 (3.5)	101 (1.8)	99.7 (3.2)	95.3 (1.3)
Ergometrinine	101 (4.2)	99.1 (1.9)	94.3 (0.7)	93.2 (4.3)	95.5 (1.7)	89.1 (2.2)	90.6 (3.9)	90.1 (4.4)	89.7 (1.9)	100 (3.5)	98.5 (1.9)	91.1 (1.9)
Ergosine	108 (2.6)	106 (5.6)	101 (3.2)	90.8 (2.0)	91.8 (2.2)	89.2 (2.6)	91.7 (2.2)	90.4 (3.1)	90.3 (1.5)	99.9 (2.7)	99.1 (3.0)	98.2 (1.1)
Ergosinine	111 (1.8)	109 (1.9)	103 (1.1)	100 (1.1)	102 (2.0)	97.7 (2.2)	92.7 (1.4)	93.6 (2.5)	93.8 (0.9)	99.2 (2.8)	98.4 (2.8)	97.5 (1.0)

Ergotamine	109 (1.9)	108 (1.7)	102 (2.8)	91.0 (2.8)	92.6 (2.8)	89.8 (3.6)	91.1 (2.2)	90.6 (3.7)	90.7 (1.3)	101 (2.9)	100 (3.1)	96.4 (2.2)
Ergotaminine	109 (1.9)	109 (0.7)	101 (0.6)	98.2 (2.0)	101 (1.5)	96.6 (1.3)	93.6 (3.5)	94.7 (1.7)	94.5 (0.6)	101 (2.3)	99.7 (1.3)	97.1 (1.5)

Table 2: Accuracy and precision analysis of fortified food samples.

CONCLUSION

This study established a workflow for the simultaneous determination of *Alternaria* toxins, ergot alkaloid epimers, and other major mycotoxins produced by the fungal genus of *Aspergillus*, *Fusarium*, and *Penicillium*. This approach can be applied to quantify multiple mycotoxins in a wide range of food commodities. A matrix-matched calibration was implemented to compensate for inherent matrix effects. The majority of analytes could be measured at a concentration as low as 0.4 µg/kg, which was satisfactory to meet the regulatory requirements.

REFERENCES

1. Commission Regulation (EU) 2015/1940 of 28 October 2015 amending regulation (EC) No 1881/2006 as regards maximum levels of ergot sclerotia in certain unprocessed cereals and the provisions on monitoring and reporting. *Off J Eur Union*. 2015;283:3-8.
2. European Commission. Summary report of the standing committee on plants, animals, food, and feed section novel food and toxicology safety of the food chain. 2021.
3. Food chemistry institute of the association of the German confectionery industry. *Alternaria* toxins: Occurrence, toxicity, analytical methods, maximum levels. 2020.
4. Naude TW, Botha CJ, Vorster JH, Roux C, van der Linde EJ, van der Walt SI, et al. *Claviceps cyperi*, a new cause of severe ergotism in dairy cattle consuming maize silage and teff hay contaminated with ergotised *Cyperus esculentus* (nut sedge) on the Highveld of South Africa. *Onderstepoort J Vet Res*. 2005;72(1):23-37.
5. Logrieco A, Moretti A, Solfrizzo M. *Alternaria* toxins and plant diseases: An overview of origin, occurrence and risks. *World Mycotoxin J*. 2009;2(2):129-140.
6. Tölgyesi Á, Kozma L, K. Sharma V. Determination of *Alternaria* toxins in sunflower oil by liquid chromatography isotope dilution tandem mass spectrometry. *Molecules*. 2020;25(7):1685.
7. Noser J, Schneider P, Rother M, Schmutz H. Determination of six *Alternaria* toxins with UPLC-MS/MS and their occurrence in tomatoes and tomato products from the Swiss market. *Mycotoxin Res*. 2011;27(4):265-271.
8. Di Mavungu JD, Malysheva SV, Sanders M, Larionova D, Robbens J, Dubrue P, et al. Development and validation of a new LC-MS/MS method for the simultaneous determination of six major ergot alkaloids and their corresponding epimers. Application to some food and feed commodities. *Food Chemistry*. 2012;135(1):292-303.
9. Krska R, Stubbings G, Macarthur R, Crews C. Simultaneous determination of six major ergot alkaloids and their epimers in cereals and foodstuffs by LC-MS-MS. *Anal Bioanal Chem*. 2008;391:563-576.
10. EFSA Panel on Contaminants in the Food Chain (CONTAM). Scientific opinion on the risks for animal and public health related to the presence of *Alternaria* toxins in feed and food. *EFSA J*. 2011;9(10):2407.
11. Commission Regulation (EU) 2021/1399 of August 2021 amending regulation (EC) No 1881/2006 as regards maximum levels of ergot sclerotia and ergot alkaloids in certain foodstuffs. *Off J Eur Union*. 2021;301:1-5.
12. Liang SH, York JL, Konschnik JD, Majer H, Steimling JA. Simultaneous determination of *Alternaria* toxins, ergot alkaloid epimers, and other major mycotoxins in various food matrixes by LC-MS/MS. *J AOAC Int*. 2023;106(2):333-340.